

Here's how to save the data on the UV-VIS spectrometers as CSV files to be opened in Excel later.

1. Select which spectrum you want to save. (The selected spectrum will have little diamonds along the line. You can only do 1 at a time.)
2. Go to "File"
3. Select "Export Selected Spectrum"
4. Select "CSV"
5. Select where you want the file to be saved. (To save it to a USB, use the drop-down menu on the bottom right of the screen)
6. Name your file, making sure it has the ".CSV" ending.
7. Select "OK".
8. Wait until the "Exporting data..." message on the bottom of the window goes away. This means your data is done exporting.